

Thermocult Incubator

Thermocult Bakterien-Brutschrank

732590

How can you regulate the temperature?

For temperature control, remove the cover cap above the main switch. Please change the temperature only with very small movements (movements clockwise increase and movements counterclockwise decrease the temperature), if necessary or desired. The temperature should not change on its own, otherwise usually the thermostat is defective.

Wie kann die Temperatur geregelt werden?

Für die Temperaturregelung die Abdeckkappe über dem Hauptschalter entfernen. Bitte die Temperatur nur mit ganz kleinen Bewegungen ändern (im Uhrzeigersinn wird die Temperatur erhöht, gegen den Uhrzeigersinn verringert), falls erforderlich oder gewünscht. Die Temperatur sollte sich nicht selbstständig ändern, andernfalls ist meist der Thermostat defekt.

EICKEMEYER®

GERMANY | EltstraÙe 8 | 78532 Tuttlingen | Deutschland | T +49 7461 96 580 0 | F +49 7461 96 580 90 | info@eickemeyer.de | www.eickemeyer.de

SWITZERLAND | Sandgrube 29 | 9050 Appenzell | Schweiz | T 071 788 23 13 | F 071 788 23 14 | info@eickemeyer.ch | www.eickemeyer.ch

UNITED KINGDOM | 3 Windmill Business Village | Brooklands Close | Sunbury-on-Thames | Surrey | TW16 7DY | UK | T 020 8891 2007 | www.eickemeyer.co.uk

DENMARK | Solbakken 26 | Hammelev | 6500 Vojens | Danmark | T +45 70 20 50 19 | info@eickemeyer.dk | www.eickemeyer.dk

NETHERLANDS | Bellweg 44 | 4104 BJ Culemborg | Nederland | T +31 345 58 9400 | info@eickemeyer.nl | www.eickemeyer.nl

ITALY | Via G. Verdi 8 | 65015 Montesilvano (PE) | Italia | T 085 935 4078 | F 085 935 9471 | info@eickemeyer.it | www.eickemeyer.it

POLAND | Al. Jana Pawła II 27 | 00-867 Warszawa | Polska | T 22 185 55 76 | F 22 185 59 40 | info@eickemeyer.pl | www.eickemeyer.pl

CANADA | 617 Douro Street, Suite #205 | Stratford | Ont. | Canada | N5A 0B5 | T +1 519 273 5558 | F +1 519 271 7114 | info@eickemeyervet.ca | www.eickemeyercanada.ca